



Connecting Devices to ProScope Mobile

Download App

- Go to the App Store and download **AirMicro** (iPod) or **AirMicro pad** (iPad)
- Turn on ProScope (power switch at base)

Network Settings

- Open the **Settings** app on all iPads you wish to connect to the ProScope
- Select **Wi-Fi**, then pick the **AirMicro** Network
(write down the number that appears after AirMicro, for example AirMicro089)
- Tap the blue arrow next to **AirMicro**
- Tap on the **Static** tab:
- In IP address: type in **192.168.X.Y**
 - The third set of numbers (X) is the number from the network name, without the zero)
 - The fourth set of numbers (Y) can be anything from 2 to 254. If multiple devices will be connecting to this ProScope, each one needs to have a different number in the Y spot.
 - For example, with a network named AirMicro098, the IP addresses for each device would be 192.168.89.2, 192.168.89.3, 192.168.89.4, 192.168.89.5 and so on.
- In Subnet Mask: type in 255.255.255.0

Open the AirMicro App

- You should be connected.
- Turn the power off and on again if you do not see the microscope.
- Be sure to turn the light on to see clearly.

When finished, be sure to switch back to your original wireless network.